



Comparative IMCStip-Based Protein A Purification on the Tecan Fluent: Impact of Resin and Elution pH on Antibody Aggregation

Todd Mullis¹ • Patrick Kates¹ • Gracie Anderson^{1,2} • L. Andrew Lee¹

¹Integrated Micro-Chromatography Systems, Inc., Irmo, SC • ²University of South Carolina, Columbia, SC

Contact: inquiries@imcstips.com

INTRODUCTION

Optimization of protein purification is critical for obtaining high recovery, purity, and monomer percentage. The optimization process is generally a time consuming and manual process. Several variables need to be accounted for including:

- Resin type
- Resin binding capacity
- Wash/Elution conditions

In this study, IMCStips are used for the purification of two antibody samples. For the optimization of protein purification, two commercially available resins were used:

- MabSelect™ Prisma: Industry standard, high binding capacity, specificity
- Praesto™ Jetted A50 HipH: Known for higher pH elution

Use of Tecan IMCStips in a fully automated antibody purification and buffer exchange process on the Tecan Fluent platform

Key findings:

- pH effect: Similar recovery at pH 3.6; higher recoveries with Praesto™ Jetted A50 HipH at pH 3.9-4.5 vs MabSelect™ Prisma
- Monomer percentage: Resin- and pH-specific effects

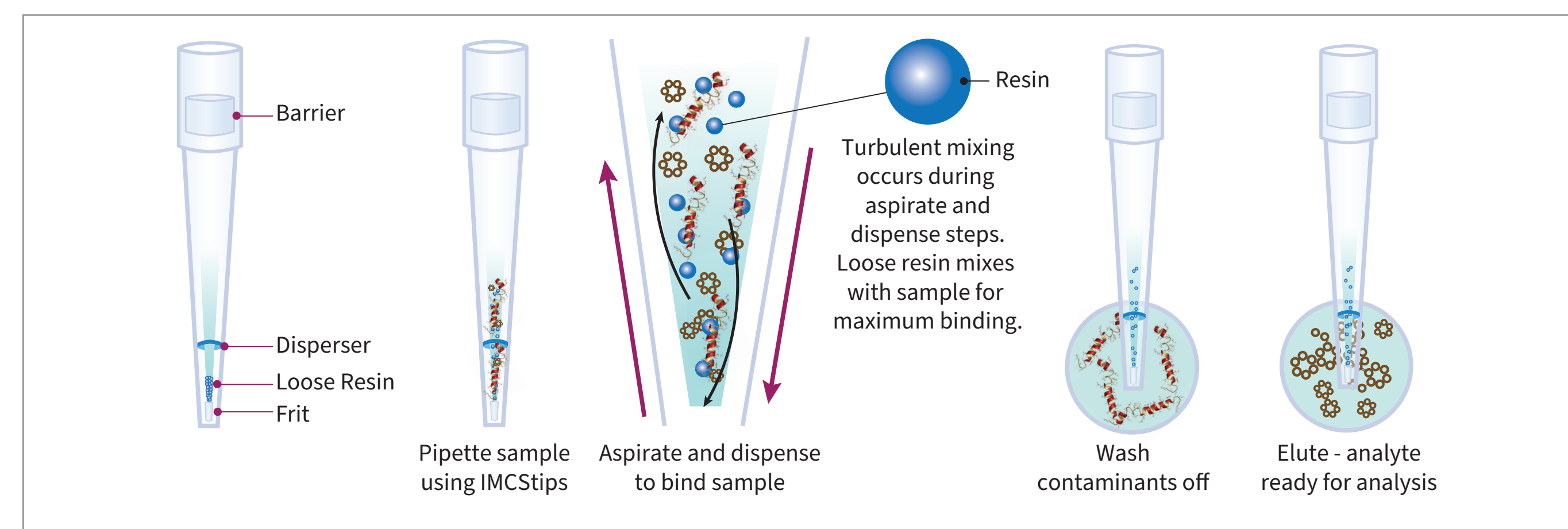


Figure 1: IMCStips® containing loose resin employ dSPE to perform efficient automated extractions.

METHODS

Fully automated Protein A purification and buffer exchange performed on the Tecan Fluent system

- 1 mL Protein A Tecan IMCStip
 - 25 µL MabSelect™ PrismaA
 - 25 µL Praesto™ Jetted A50 HipH
- 1 mL SizeX₁₅₀ Tecan IMCStip

Monoclonal antibodies captured from clarified cell culture supernatants using MabSelect™ PrismaA and Praesto™ Jetted A50 HipH

Elution pH range: 2.9 to 4.9

Analyzed purified antibody samples for recovery and monomer percentage using:

- Absorbance at 280 nm
- SDS-PAGE
- Size-exclusion chromatography (SEC)

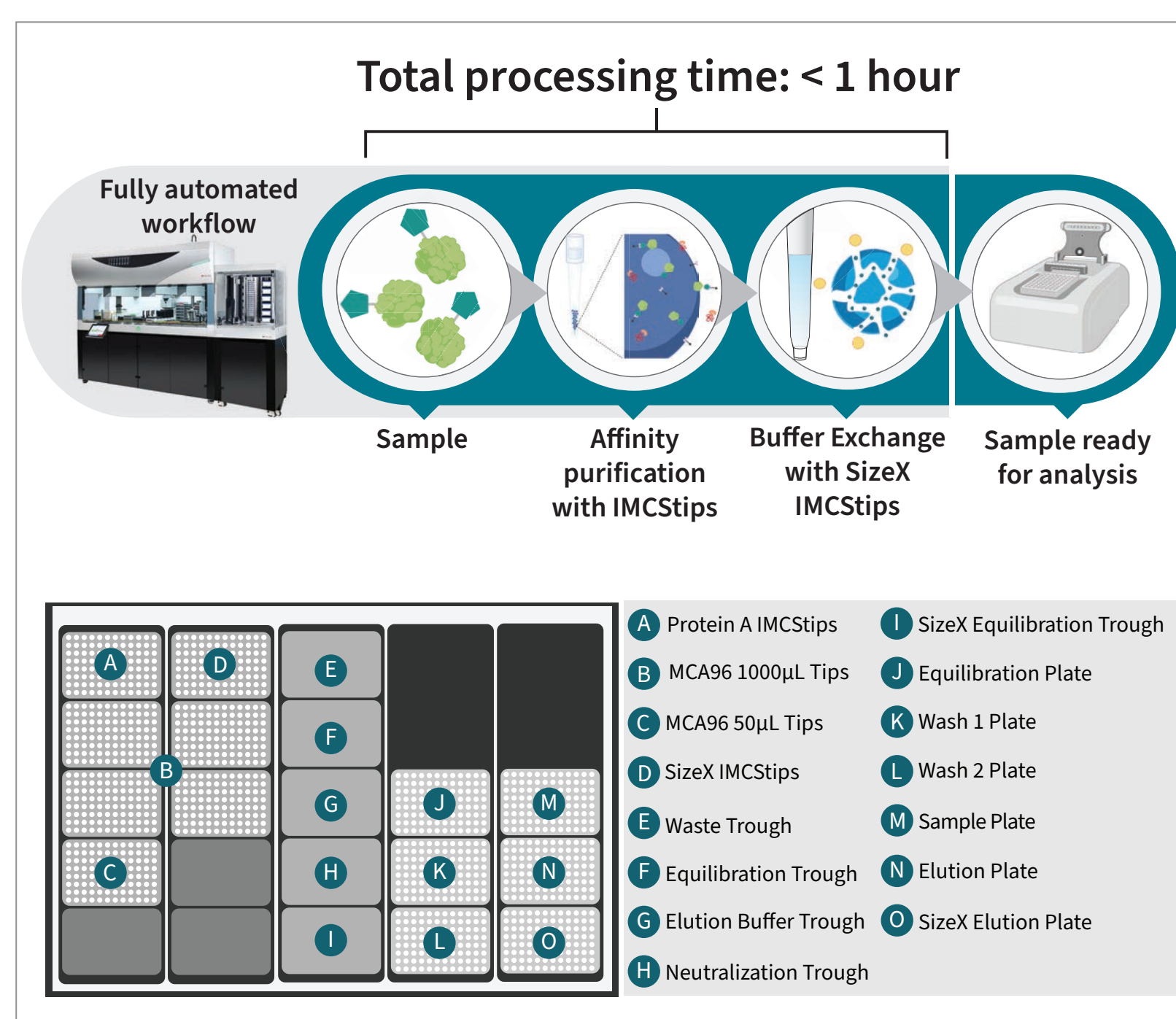


Figure 2: Sample deck layout for automated protein purification followed by buffer exchange of 96 samples using IMCStips on the Tecan Fluent.

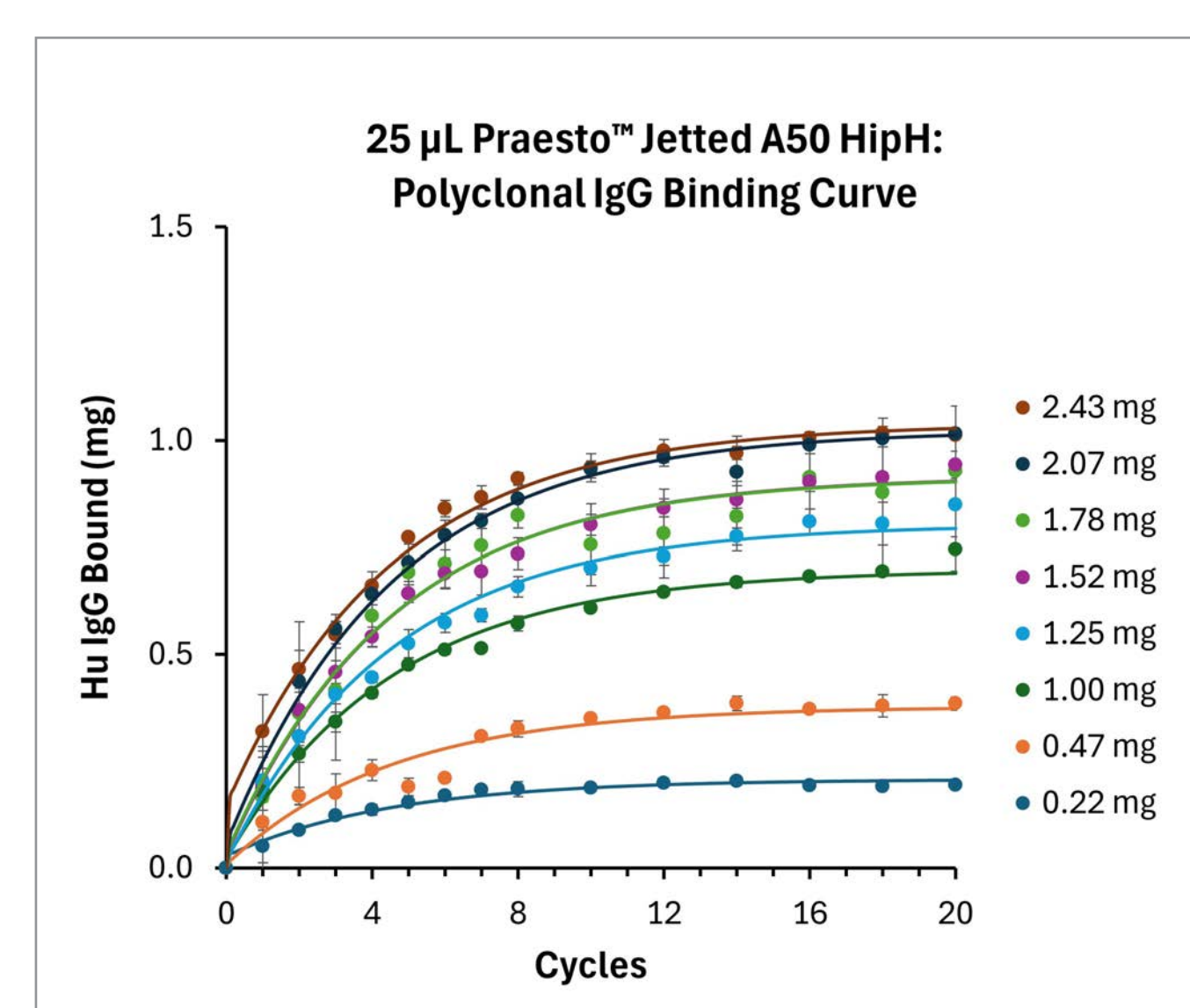


Figure 3: Binding kinetics for 25 µL HipH 1 mL Tecan IMCStips.

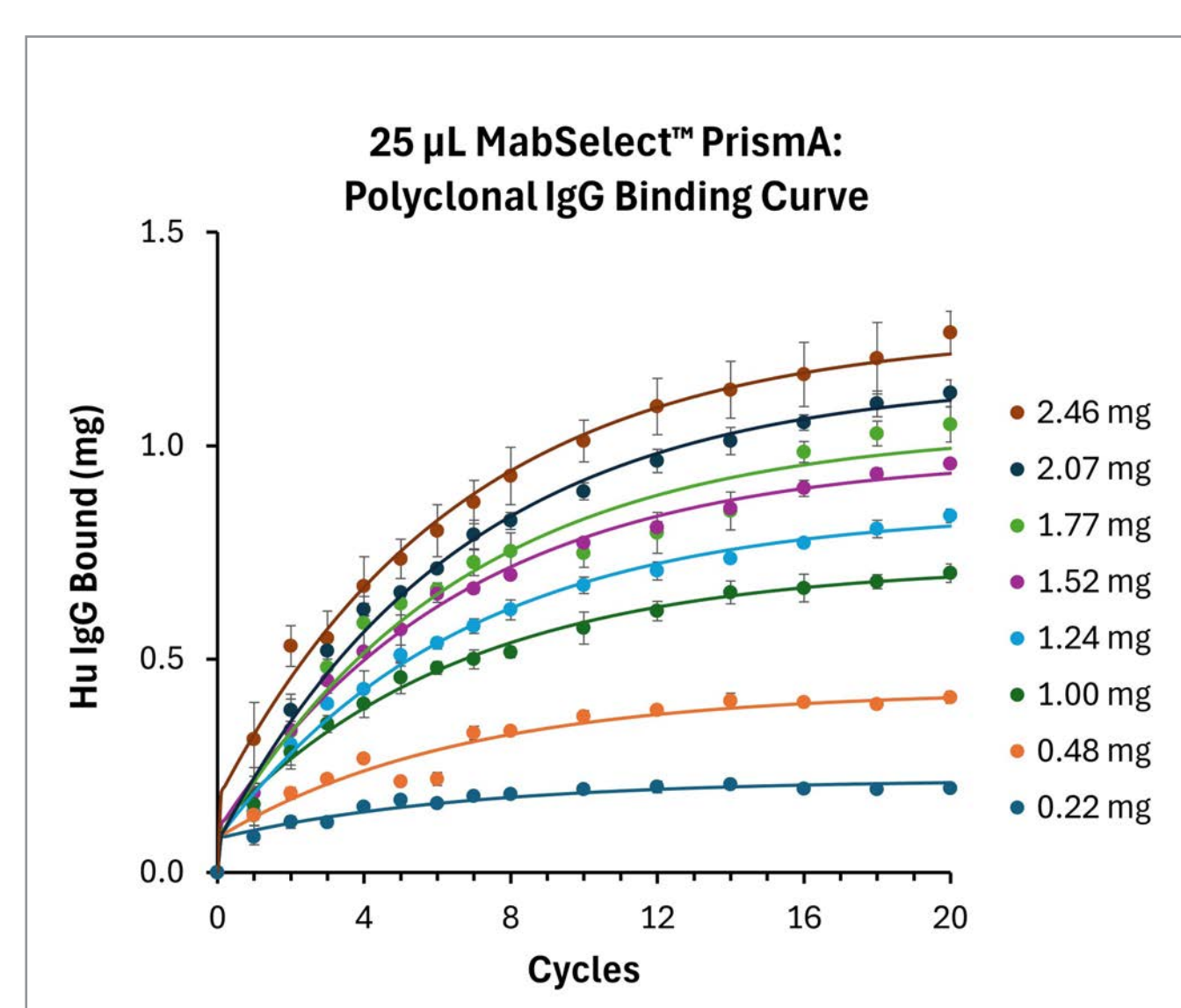


Figure 4: Binding kinetics for 25 µL Prisma 1 mL Tecan IMCStips.

MabSelect™ Prisma v Praesto™ Jetted A50 HipH: Recovery Comparison

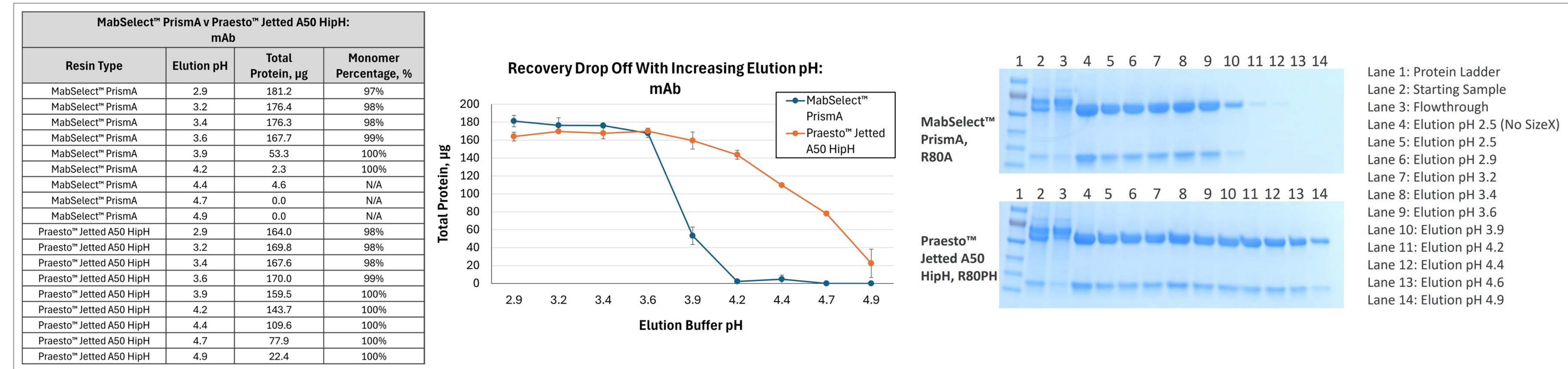


Figure 5: Comparison of recovery between MabSelect™ PrismaA and Praesto™ Jetted A50 HipH with different elution buffer pH's. For the purification, a monoclonal antibody expressed in cell culture supernatant was used.

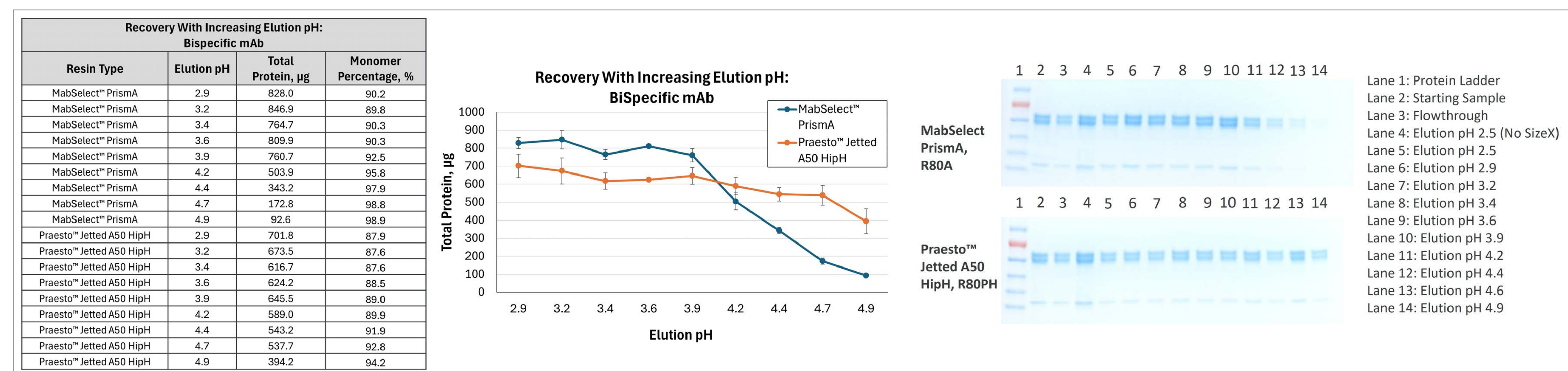


Figure 6: Comparison of recovery between MabSelect™ PrismaA and Praesto™ Jetted A50 HipH with different elution buffer pH's. For the purification, a bispecific monoclonal antibody expressed in cell culture supernatant was used.

mAb: MabSelect™ Prisma

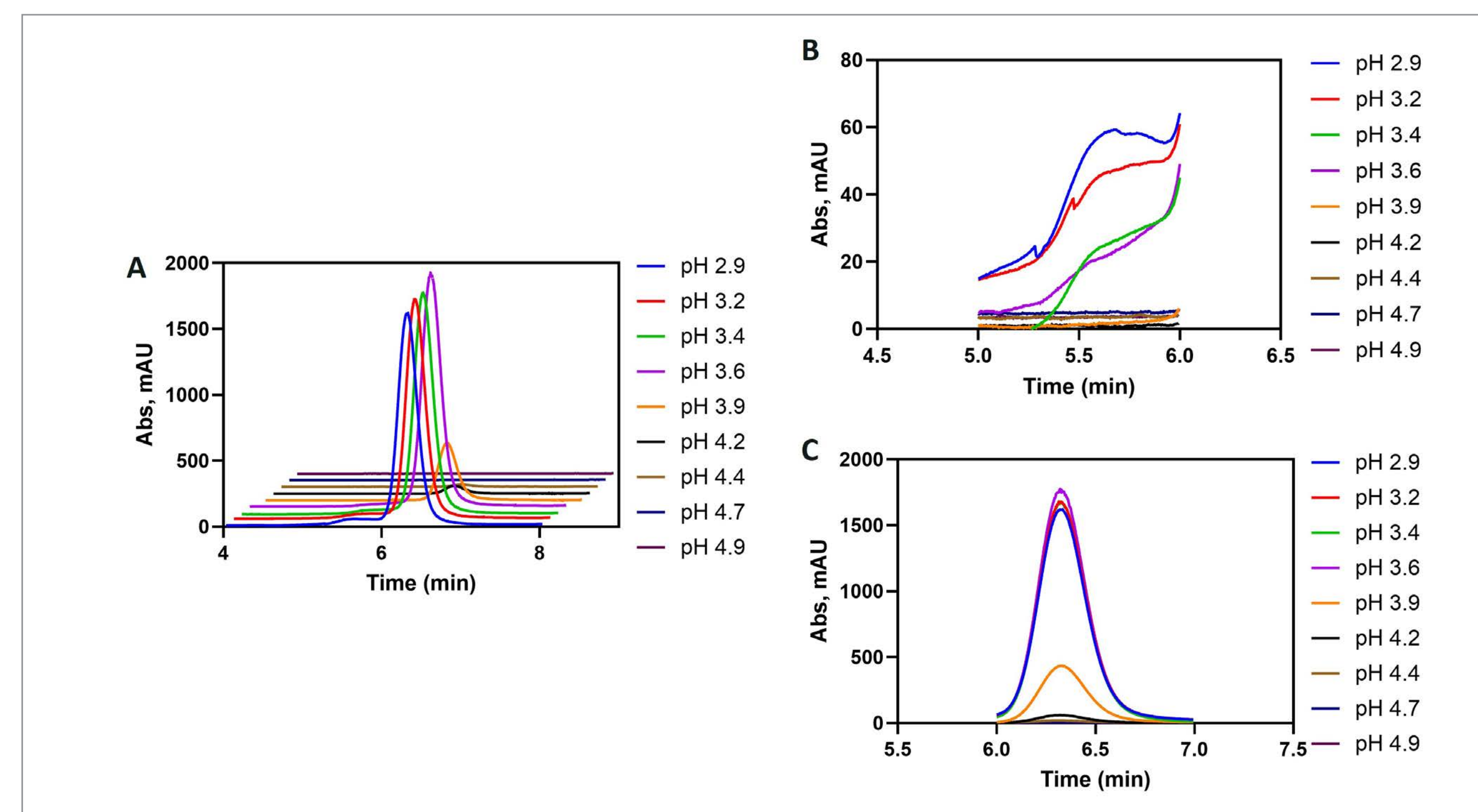


Figure 7: Comparison of chromatography from MabSelect™ Prisma purified mAb eluted at different pHs. A) Full chromatograph, B) Aggregate, C) Monomer.

mAb: Praesto™ Jetted A50 HipH

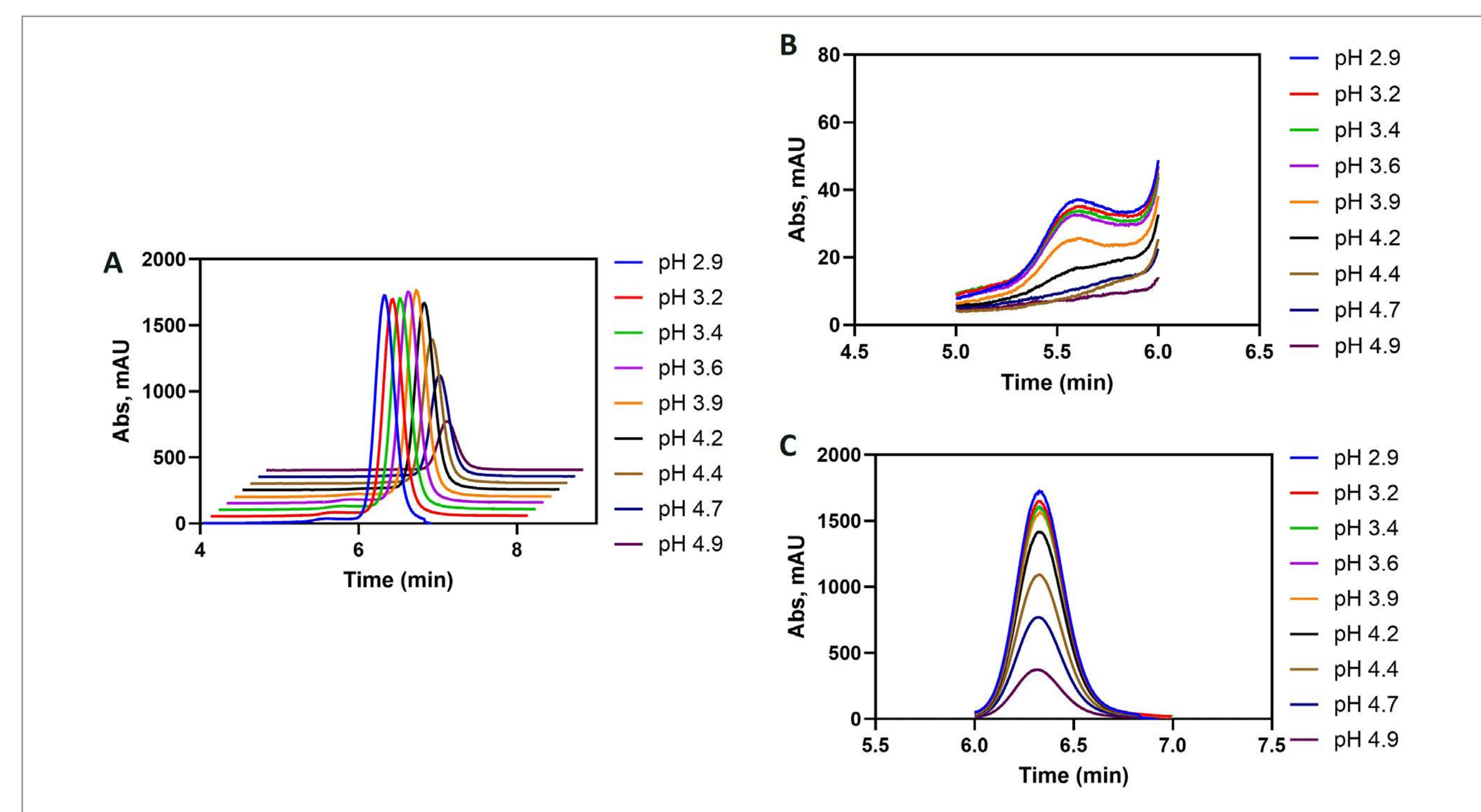


Figure 8: Comparison of chromatography from Praesto™ Jetted A50 HipH purified mAb eluted at different pHs. A) Full chromatograph, B) Aggregate, C) Monomer.

RESULTS

- Automated method established on the Tecan Fluent enables purification and buffer exchange of 96 samples in under 60 minutes.
- MabSelect™ Prisma has a higher binding capacity than Praesto™ Jetted A50 HipH.
- Using two different cell supernatant samples containing mAbs:
 - Both resins achieve similar recovery across a range of low pH values.
 - Comparable yields up to pH 3.6.
- At low pH elution conditions:
 - Monomer percentages of the two antibodies tested were notably influenced by the resin used even at the same elution pH.
 - As pH increases, monomer percentages increases.
- At higher pH elution conditions (3.9 and above):
 - Praesto™ Jetted A50 HipH resin yields significantly higher recovery.

CONCLUSION

- IMCStips allow for efficient optimization of protein purification allowing for streamlined testing of resin types, binding capacity, and elution conditions.
 - Traditional methods such as FPLC would require significantly more hands-on time.
- Data shows that recovery and monomer content are influenced by:
 - The specific resin used for purification.
 - The elution pH.
- This work paves the way for automating and optimizing antibody purification screening strategies to improve both yield and monomer content.

DOWNLOAD POSTER

Scan QR code to get access to this poster online.

